

Intrinsic Dissolution Apparatus

Operator's Manual

Limitation of Liability

The information in this document is subject to change without notice. Dissolution Accessories makes no warranty of any kind with regard to this material, including, but not limited to, the implied warranties or merchantability and fitness for a particular purpose. Dissolution Accessories shall not be liable for errors contained herein or for incidental consequential damages in connection with the furnishing, performance, or use of this material. All rights are reserved. Reproduction, adaptation, or translation without prior written permission is prohibited, except as allowed under copyright laws.

Table of Contents

1. Introduction	4
History of the Intrinsic Dissolution Apparatus	4
2. Setup	6
Unpacking Your Intrinsic Dissolution Apparatus	6
Assembling the Intrinsic Apparatus	8
3. Operation	12
Sample Preparation and Compaction Procedure	12
Suggested Equipment	12
Suggested Materials	12
Preparation of the Compact	13
General Test Procedure	14
Sampling and Analysis	15
Calculation and Interpretation	15
4. Maintenance	16
Cleaning	16
Index	17

List of Figures

FIG. 1. Intrinsic Dissolution Apparatus	5
FIG. 2. Intrinsic Dissolution Apparatus Items	6
FIG. 3. Assembly of the Intrinsic Apparatus	7
FIG. 4. Upside Down View of Surface Plate with Recessed Screws	8
FIG. 5. Center Opening of Die	8
FIG. 6. Inserting the Punch in the Intrinsic Apparatus	7
FIG. 7. Tighten Slotted Nylon Screws	10
FIG. 8. Holder Attached to the Die	10
FIG. 9. Intrinsic Apparatus Attached to the Dissolution Tester	11
FIG.10. Experimental Data Plot	15

1. Introduction

The measurement of intrinsic dissolution rates is an increasingly important tool in the Pharmaceutical research and development laboratory. It allows the characterization of both the crystallized forms and polymorphs of pure drugs and drug formulations by exposing a constant surface area to the dissolution medium. When measuring intrinsic dissolution rates, the parameters which control the rate of dissolution, such as surface area exposed to the medium, temperature, stirring rate, and pH, are kept constant.

Additional information is available in the USP General Chapters, Intrinsic Dissolution <1087>.

This product is warranted to be free from manufacturing defects at the time of delivery. Please report any problems upon receipt of merchandise to info@dissolutionaccessories.com

History of the Intrinsic Dissolution Apparatus

The evolution of pharmaceutical testing techniques led to the development of what has become known as the Wood Apparatus for intrinsic dissolution testing by Dr. John Wood in 1963. The original Wood Apparatus consists of a punch and a die fabricated of type 316 stainless steel. The base of the die has three threaded holes for the attachment of a polished steel plate which provides a mirror-smooth base for the compaction of the drug substance under study. To determine a material's intrinsic dissolution rate, a measured amount of the material is placed into the 0.8 cm diameter cavity of the die. The punch is inserted into the die cavity on top of the test material. The material is compressed using a benchtop tablet press (a hole in the head of the punch allows insertion of a metal bar to facilitate removal after the test). A pellet of the compressed material, called a compact, is formed in the cavity with a single face of defined area (0.5 cm²) exposed on the bottom of the die. The bottom of the die cavity is threaded so that at least 50% of the compact can dissolve without falling out. The top of the die has a threaded shoulder which allows it to be attached to a holder. The holder is mounted on a laboratory stirring device, and the entire die with the compact still embedded is immersed in the dissolution medium (often contained in a 400 mL water-jacketed beaker) and rotated by the stirring device. Samples are taken and assayed at periodic intervals, enabling the calculation of the dissolution rate per centimeter squared.

In our design, the walls of the die cavity are grooved for holding the compressed drug substance within the die cavity during the test. The ends of the cavity are perfectly squared off and the surface area exposed remains constant as the compact dissolves. High pressures are sometimes required to form the compact and the type 316 stainless steel originally used could deform. The Varian design uses hardened steel, which stands up to high pressure, and all surfaces in contact with the drug substance are polished for even compression. The bottom of the die contains three threaded holes, enabling the surface plate to be screwed onto the die to prevent slippage during pressing.

Due to the weight of the intrinsic device, the Intrinsic Dissolution Apparatus is designed for use with a dissolution tester with a secure shaft mounting system such as a chuck and collet. The dissolution apparatus offers more precise speed and water bath temperature control than a benchtop stirrer and water-jacketed beaker. Also, multiple tests can be run simultaneously.

Note

The punch is made from hardened steel. Both the punch and the surface plate should be oiled when not in use. See Chapter 5, "Maintenance," on page 27. All other parts on the intrinsic dissolution device are fabricated from type 316 stainless steel.

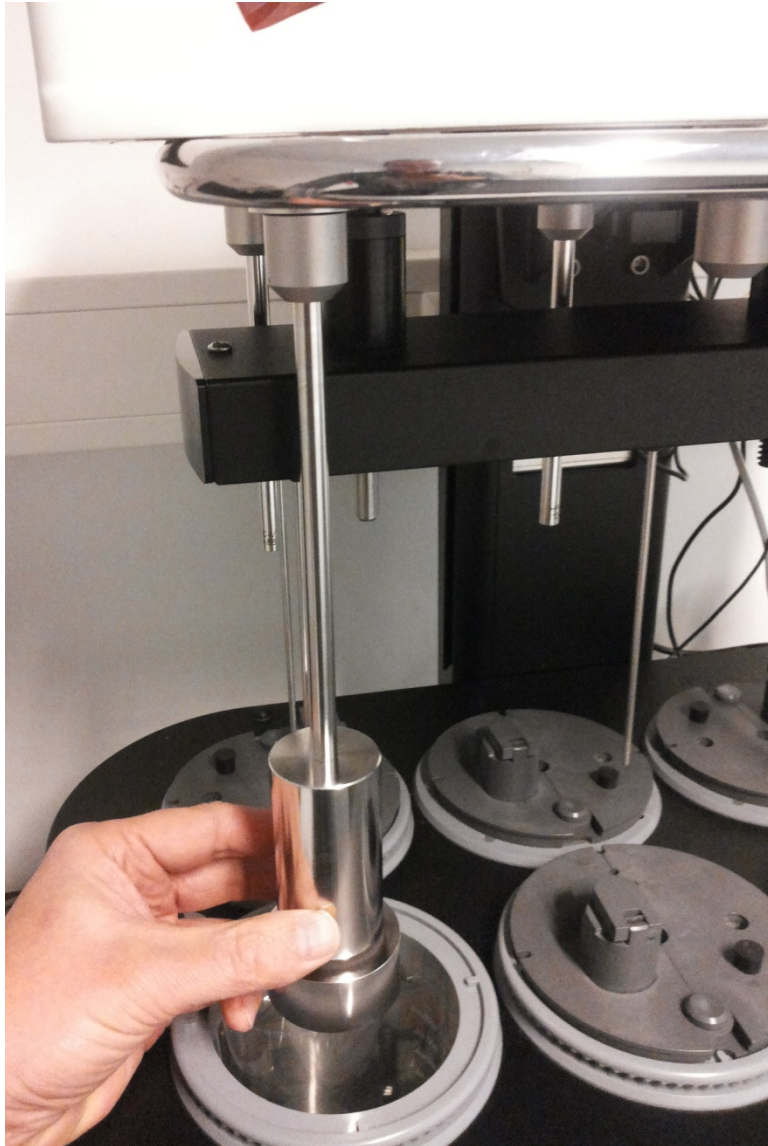


FIG. 1. Intrinsic Dissolution Apparatus

2. Setup

Unpacking Your Intrinsic Dissolution Apparatus

- Step 1 Carefully remove the apparatus from its shipping carton. The items are heavy; hold them firmly to prevent them from dropping.
- Step 2 Be sure to remove all parts before discarding or storing the packaging.
- Step 3 Place the items on a clean, dry, level section of the bench top or table.



FIG. 2. Intrinsic Dissolution Apparatus Items

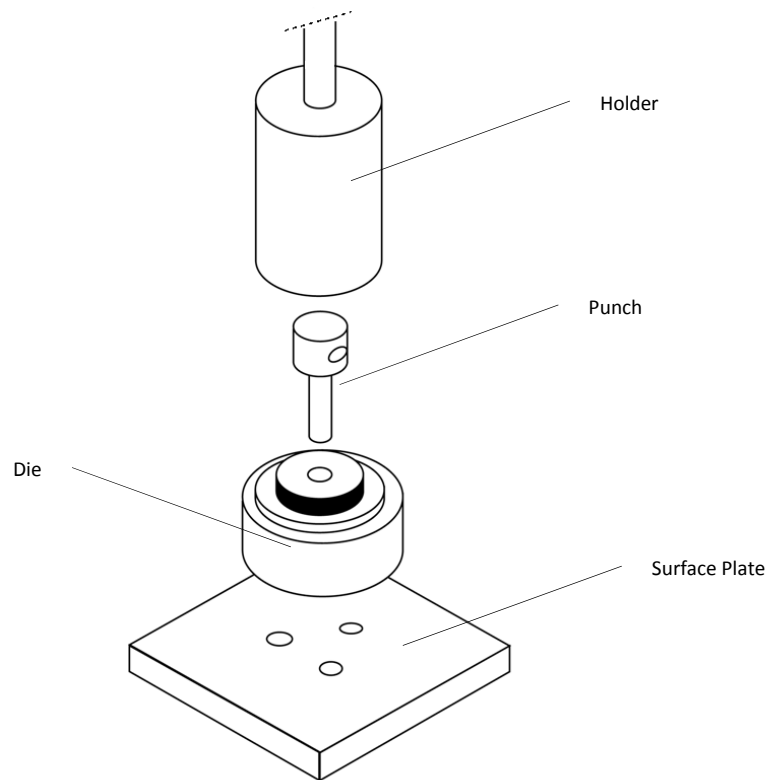


FIG. 3. Assembly of the Intrinsic Apparatus

Note

The punch and surface plate are made from hardened steel. Both the punch and the surface plate should be oiled when not in use. See chapter 4, "Maintenance," on page 14. All other parts on the intrinsic dissolution apparatus are fabricated from type 316 stainless steel.

Assembling the Intrinsic Apparatus

- Step 1 Ensure all traces of oil are removed from the punch and surface plate.
- Step 2 Place the surface plate on a flat surface so that the recessed holes are on the bottom.

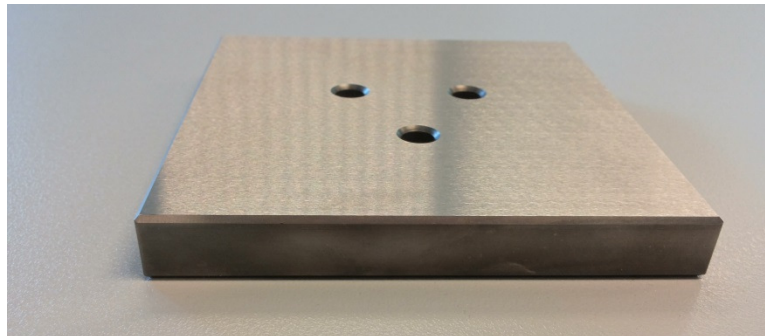


FIG. 4. Upside Down View of Surface Plate with Recessed Screws

- Step 3 Place the die on top of the surface plate aligning the three holes in the die with the three holes in the surface plate.
- Step 4 Pick up the assembly holding the two pieces in the appropriate alignment. Turn the assembly over and use the fastening screws to secure the pieces together.
- Step 5 Use an Allen key to tighten the fastening screws. Turn the assembly over again and place it on a laboratory table. Because the holes on the bottom of the surface plate are recessed, the surface plate sits flat when the screws are tightened.
- Step 6 Place the washer on the die.
- Step 7 Accurately weigh out the test compound.
- Step 8 Carefully transfer the test compound into the cavity of the die.

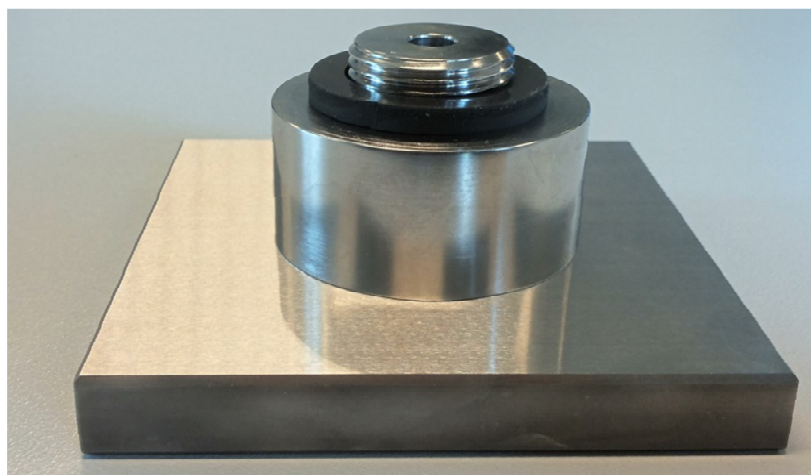


FIG. 5. Center Opening of Die

Step 9 Place the punch in the cavity over the test compound.

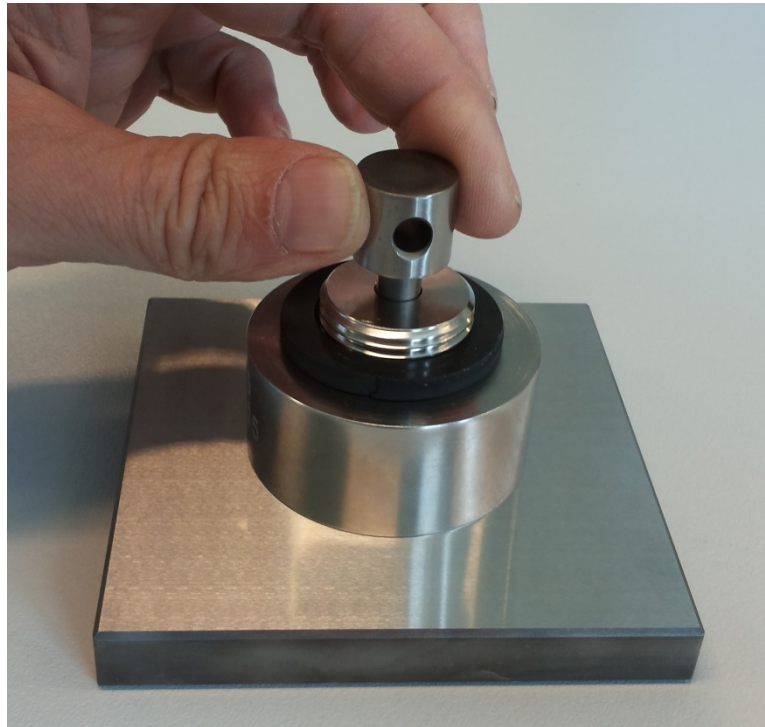


FIG. 6. Inserting the Punch in the Intrinsic Apparatus

- Step 10 Place the entire assembly in a laboratory press equipped with an accurate pressure gauge.
- Step 11 Apply pressure to compress the material. Take care not to over compress the drug substance.
- Step 12 Remove the assembly from the laboratory press.
- Step 13 Turn the assembly over and remove the three fastening screws from the surface plate. Be careful not to let the punch and die fall when the fastening screws are removed.
- Step 14 Remove the punch and die from the surface plate.
- Step 15 Place the slotted nylon screws provided with the apparatus in the holes in the bottom of the die. Use a screwdriver to tighten the screws.

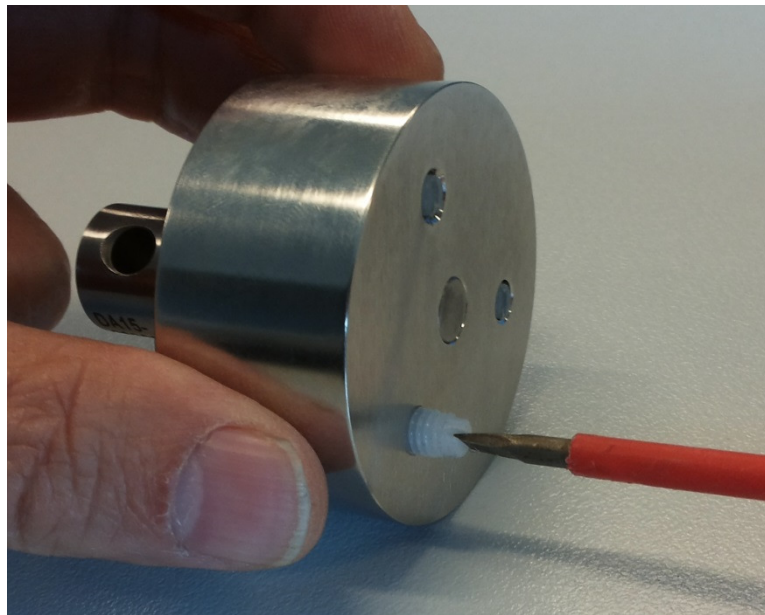


FIG. 7. Tighten Slotted Nylon Screws

Step 16 Insert the top of the die, with the punch still in place, into the holder and tighten.

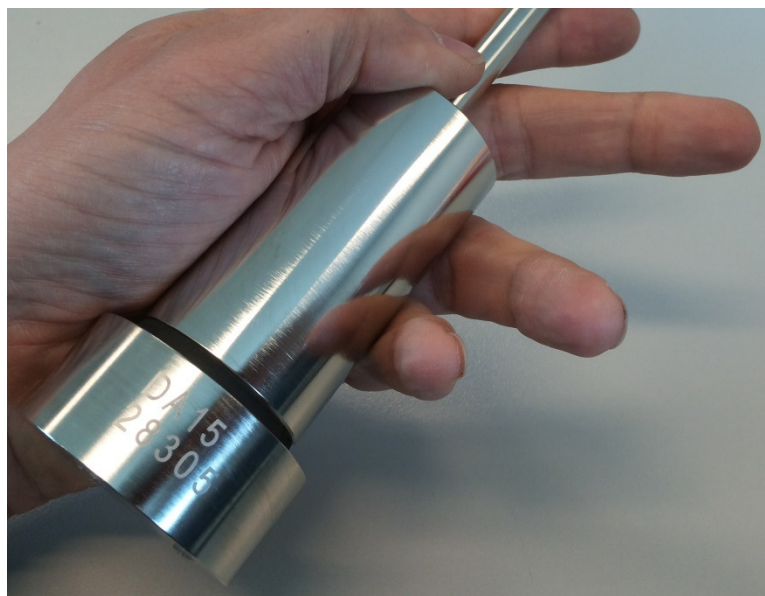


FIG. 8. Holder Attached to the Die

Step 17 Remove all loose powder from the surface of the die.

Step 18 Insert the intrinsic assembly into the dissolution apparatus spindle assembly and tighten. Carefully lower the drive unit and adjust the height so that the exposed surface of the compacted pellet is 3.8 cm from the bottom of the vessel. See Fig. 9, "Intrinsic Apparatus Attached to the Dissolution Tester," below.

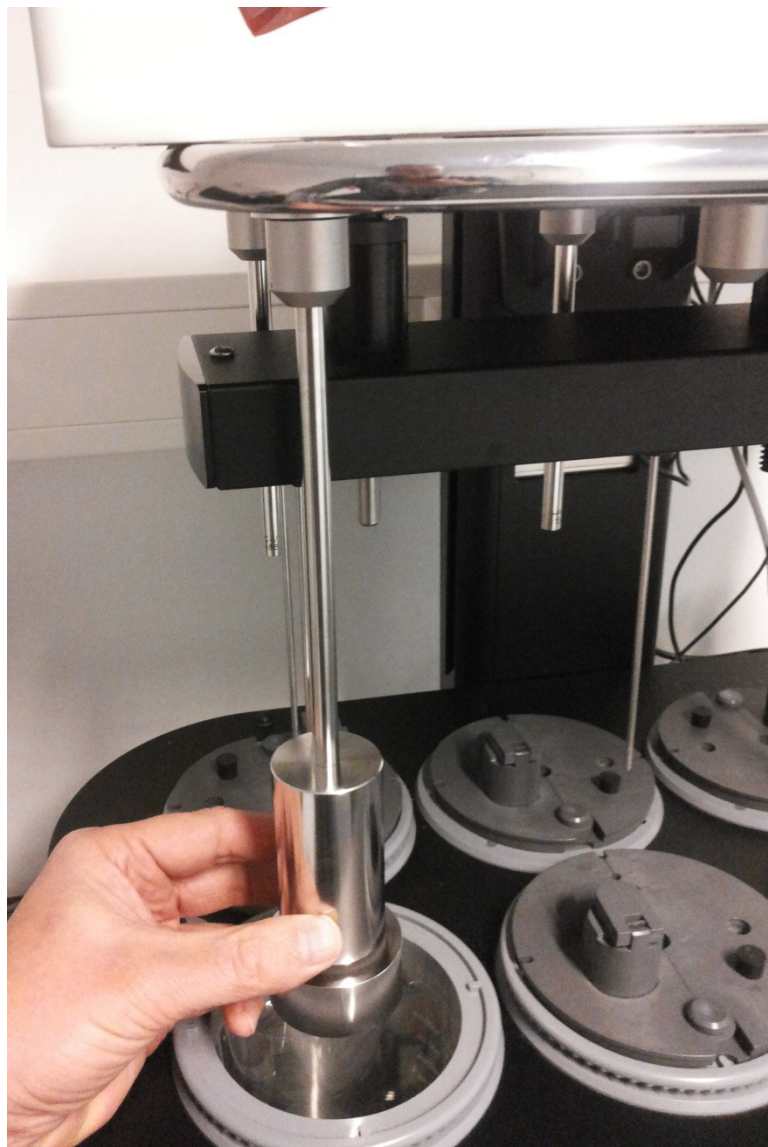


FIG. 9. Intrinsic Apparatus Attached to the Dissolution Tester

Note

The dissolution apparatus should have suitable means of securing the intrinsic device. Due to the weight of the device, a chuck and collet system works best.

3. Operation

Sample Preparation and Compaction Procedure

Note

The following procedures are intended for illustration and guidance only. Actual practice depends upon the nature of the material being tested and the equipment used.

Suggested Equipment

- Intrinsic Dissolution Assembly (including punch, die, holder and surface plate)
- Hydraulic bench top press with accurate pressure gauge (Carver Press or E-Z Press)
- Dissolution tester with variable speed control
- Amber (tinted) vessels for light-sensitive compounds
- Sampling device (syringe with cannula or sampler)
- Detection device (UV-Vis spectrophotometer or HPLC)

Suggested Materials

- Test compound
- Degassed dissolution medium
- Standard solution prepared with same degassed dissolution medium

Preparation of the Compact

- Step 1 Accurately weigh out the compound to be tested on a piece of tared weighing paper.
- Step 2 Attach the surface plate to the underside of the die and secure with the three fastening screws provided.
- Step 3 Carefully and evenly pour the weighed material into the die cavity.
- Step 4 Place the punch into the cavity.
- Step 5 Place the intrinsic assembly into a suitable hydraulic press.
- Step 6 Compress the material in a hydraulic press for one minute at the minimum compression pressure necessary to form a non-disintegrating compacted pellet or compact.

Caution

Use no more than 5,000 pounds of force on the die. Even though it is fabricated of hardened steel, excessive force will jam the punch into the die cavity. Damage caused by excess force is not covered under warranty.

- Step 7 Detach the surface plate and screw the die with punch still in place into the holder.
- Step 8 Remove all loose powder from the surface of the die by blowing it with compressed air or nitrogen.

General Test Procedure

- Step 1 Raise the drive unit to its highest position.
- Step 2 Loosen the hex screw of the end knob or turn chuck on one of the spindle assemblies by turning it counterclockwise. Carefully insert the shaft into the spindle assembly. Because the shaft is held very firmly, it could be necessary to use moderate force.
- Step 3 Adjust the height so the exposed surface of the compact will be 3.8 cm from the bottom of the vessel when the drive unit is lowered. Tighten the hex screw of the end knob or chuck to hold it in place.
- Step 4 Repeat the procedure for the remaining intrinsic assemblies. Leave the drive unit in the up position.
- Step 5 Set the spindle speed for the desired rate of dissolution. Speeds of 50 - 100 RPM are typical.
- Step 6 Lower the intrinsic apparatus into the vessel containing the medium at 37 °C, as determined by the Arrhenius equation, until the assembly is at the proper depth.

Note

When lowering the assembly into the media, ensure there are no air bubbles on the exposed surface of the compact. If air bubbles appear, tap the shaft gently to dislodge them, taking care not to damage the compact.

- Step 7 Start the spindle rotation at time zero.
- Step 8 Pull samples at the appropriate time intervals.
- Step 9 Upon completion of the test, remove the assembly from the medium and follow the cleaning instructions in 4, "Maintenance," on page 16.

Sampling and Analysis

Take samples at preset time intervals over the predetermined length of the test. The total test length will vary with the nature of the material. In all cases, samples should be taken until a minimum of ten percent of the test material has gone into solution. Prepare a standard curve and assay samples using a suitable means of detection.

Calculation and Interpretation

An integrated form of the modified Noyes and Whitney equation is used for calculation. As long as sink conditions are maintained, the concentration gradient can be considered a constant. More detailed information can be found in *Dissolution, Bioavailability, and Bioequivalence* by Hamed M. Abdou.

The data for the cumulative amount dissolved at each time point should be corrected for sampling losses. To calculate the intrinsic dissolution rate, plot the cumulative amount of sample dissolved per unit area of the compact against time until 10% is dissolved. The cumulative amount dissolved per unit area is given by the cumulative amount dissolved at each time point divided by the surface area exposed (0.5 cm^2). Linear regression should then be performed on data points up to and including the time point beyond which 10% is dissolved. The intrinsic dissolution rate of the test material (in $\text{mg}/\text{min}/\text{cm}^2$) at the selected stirring speed is given by the slope of the regression line.

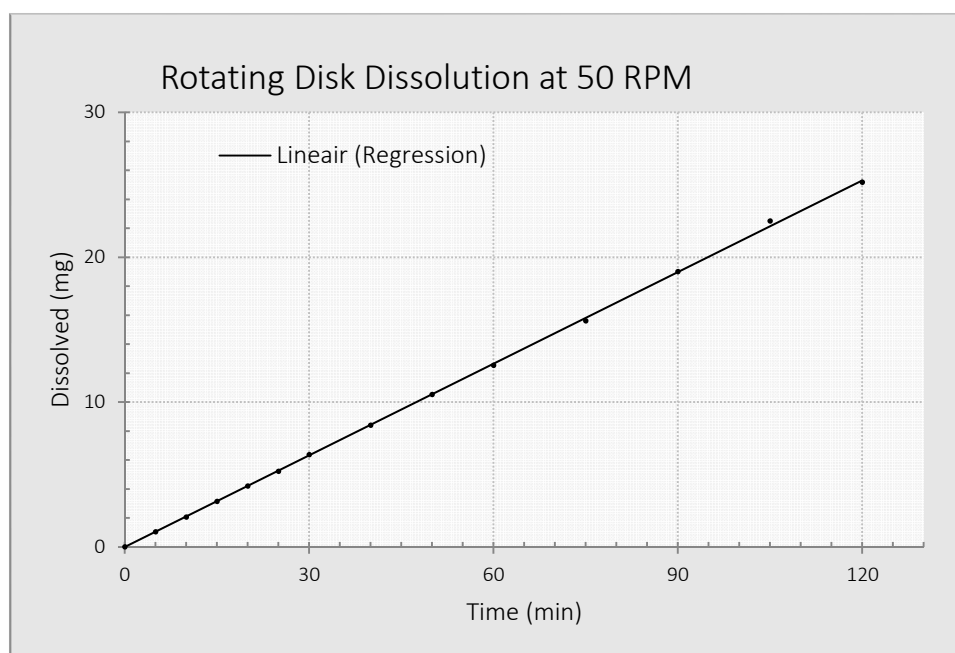


FIG. 10. Experimental Data Plot

4. Maintenance

Cleaning

After each use, it is necessary to clean the Intrinsic Dissolution Apparatus. Complete the following steps:

- Step 1 Thoroughly rinse the die, punch, holder, and surface plate with water.
- Step 2 Dry completely.
- Step 3 Apply a light coating of light machine oil to the punch and surface plate to prevent rust and corrosion.

Caution

Take careful steps to clean and lubricate your Intrinsic Dissolution Apparatus or the punch and surface plate will rust and become severely pitted.

Index

a

analysis 15
assembling the intrinsic apparatus 8

c

calculation 15
cleaning instructions 16
compaction procedure 13

d

dr. john wood 4

e

equipment, suggested 12
experimental data plot 15

g

general test procedure 14

h

history 4

i

interpretation 15
introduction 4

m

maintenance 16
materials, suggested 12

o

operating your intrinsic dissolution
apparatus 12

p

preparation of the compact 13

s

sample preparation 13
sampling 15
setting up the intrinsic dissolution apparatus 11

u

unpacking your intrinsic dissolution apparatus 11

w

wood apparatus, history 4
wood, dr. john 4